Sartorius-Wedge Quick Start

([German Version](#) p. 17)

Introduction	2
Connecting the Balance to your PC.....	2
Determining the COM Port that the Balance is Connected To	5
Modifying the Baud Rate in WinWedge	6
Inputting Balance Data into Other Windows Programs	7
Inputting Balance Readings into Any Windows Program	7
Collecting Balance Readings Using a Ready-Made Excel Workbook	8
Inputting Balance Readings into Any Excel Workbook.....	9
Connecting a balance to an Ethernet network	11
Example - Configuring the Network Settings in a Cubis Balance	12
Configuring TCPWedge to Connect to the Balance	13

TAL Technologies, Inc.

2101 Brandywine Street, Suite 102, Philadelphia PA 19130 • 215-496-0222 • www.taltech.com

© 1994 – 2015 • All Rights Reserved • Revision 1c • Last Updated 2/24/2015

Introduction

This document is designed to show how to interface your Sartorius Balance to a PC using the Sartorius-Wedge software. The Sartorius-Wedge software is a tool that allows you to input balance data directly into any Windows application program including Excel, Access, most Laboratory Information Management System (LIMS) software, etc.

Connecting the Balance to your PC

Most Sartorius balances are available with a RS232 serial port or a USB port for connecting the balance directly to a PC. Some models are also available with an Ethernet port for connecting the balance directly to a network. In the first part of this document we will show how to connect the balance directly to a PC using either the RS232 serial port or the USB port and the second section will show how to connect the balance to your network.

The RS232 serial ports on a Sartorius balance will normally have a 25-pin D-type connector and the USB port will be a USB Type-B port. The connectors are located on the back of the balance and look like the following connectors:

25-Pin RS232 Female Connector

Type B USB Port

After you determine what connectors are available on your balance, the next thing to do is to look on your PC to see what types of ports it has available. The RS232 serial ports on most PCs have 9-Pin male connectors and look like this:

Almost all PCs built within the last 15 years will have several USB ports with Type-A ports that look like the following:


Type A USB Port

If your balance has a USB port and your PC also has a USB port then this will probably be the best way to connect the two using a standard USB cable with a USB Type-A connector on one end and a USB Type-B connector on the other. When you connect the balance to your PC for the first time using a USB cable connection, Windows will attempt to install a device driver for the balance. If your PC is connected to the Internet, the driver should get installed automatically but if you are not connected to the internet, you will need to install the drivers manually using either a driver CD that came with your balance or by downloading the drivers from the following web site:


<http://www.ftdichip.com/FTDrivers.htm>

If your balance only has an RS232 serial port then this is your only option for connecting it to a PC. If your PC does not have any RS232 serial ports but it does have available USB ports then you can purchase an add-on RS232 to USB serial adapter from Sartorius. They are inexpensive and easy to install. When you first insert the adapter into your PC's USB port and the PC is connected to the Internet, Windows will automatically download and install a driver that is needed for the adapter to work. If you are not connected to the internet then you will need to install the driver manually using the driver installation CD that was provided with the adapter.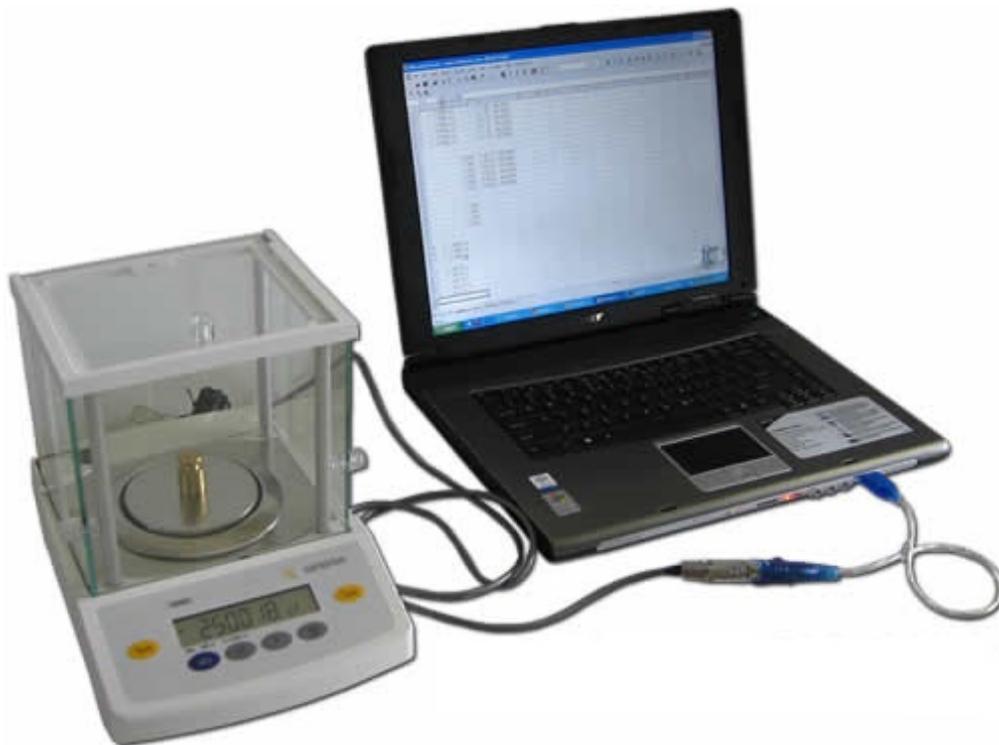

Determining the COM Port that the Balance is Connected To

Whether you connect your balance to a USB port or to a built in RS232 serial port or to a RS232 to USB adapter, as far as your PC is concerned, the balance will appear to be connected to a RS232 serial port. The next step therefore is to determine what COM port the balance is connected to. The way to do this is to open the Windows Device Manager and look in the section labeled “Ports (COM & LPT)”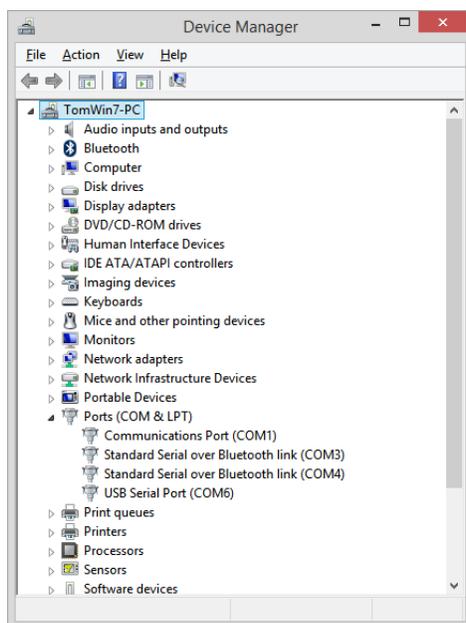

If your PC has a built in COM port, it will normally show up as either COM1 or COM2 in the Device Manager. If you connect the balance to a USB port – either directly or through a RS232 to USB adapter, the COM port will typically appear as COM3 or greater.

Once you know what COM port the balance is connected to, you can either use the Excel workbook that was provided with the Sartorius-Wedge software to input weight values into an Excel worksheet or you can manually configure the WinWedge component of the Sartorius-Wedge software to input the data from your balance into any Windows application program that you like.

In order to capture data from the balance through a serial or USB connection, you will need to know the COM port that the instrument is connected to and also the Baud Rate, Parity and Data Bits settings that the instrument is configured to use. All of the latest Sartorius instruments are shipped pre-configured with the factory default settings of 9600 Baud, Odd Parity and 8 Data Bits. Older instruments may have used 1200 Baud, Odd Parity and 7 Data Bits.

Note: It is possible to change the serial communications parameters that the balance uses. You can check to see what the settings are from the menu options available on the balance. If you think that the parameters are not set to their default settings, please consult the user’s manual for the particular instrument that you are using to learn how to navigate the menus and either check or set the serial communications parameter. The Sartorius-Wedge software is pre-configured to use the serial communications settings of 9600 Baud, Odd

Parity and 8 Data Bits therefore we recommend that you make sure that your balance is also configured to use these same settings.

The Sartorius-Wedge software is a specialized version of a program called WinWedge Professional. WinWedge is a software tool that allows you to input data from instruments connected to either a serial port or a TCP/IP network. It is primarily used to feed data from the instrument directly into some other application program that does not directly support data input through a serial or network port. In its simplest mode, WinWedge works by stuffing the data from your balance into the Windows keyboard input stream thereby making it appear as if data from the balance is actually being typed in on the keyboard. WinWedge has many powerful features that you can use to build very sophisticated data collection applications that work directly with any other Windows program. In order to use WinWedge effectively for any non-trivial application, we strongly recommend that you read the WinWedge users manual thoroughly to learn about all of its features and capabilities.

Modifying the Baud Rate in WinWedge

The Sartorius-Wedge software comes with a number of WinWedge configuration files that are used with the included sample Excel workbooks. The configuration files were created assuming that the balance you are using has been configured to use the serial communications parameters of 9600 Baud, Odd Parity and 8 Data Bits. If you are using different settings in your balance, you will either need to configure your balance to use these same settings or you will need to modify the settings in the WinWedge configuration files. The provided WinWedge configuration files are named BalanceCOMn.SW3, SARTOn.SW3 or MARK3-n.SW3. The number “n” in the file names represents the COM port that the particular configuration file has been configured for. To change the baud rate in any of these files, run WinWedge.exe and select File – Open from the WinWedge main menu and open the configuration file with the number “n” that matches the COM port that your balance is connected to. Next, you would select Port – Settings from the WinWedge main menu and change the serial communications settings in the window that appears to match the settings for your balance. Finally, you would select File – Save from the WinWedge main menu to save the new configuration settings and then File – Exit to exit WinWedge.

Inputting Balance Data into Other Windows Programs

The Sartorius-Wedge software comes with several different tools that are designed to make it quick and easy to start inputting data from a Sartorius balance directly into any Windows program as well as a few tools that are specifically designed for capturing balance data into Excel.

Inputting Balance Readings into Any Windows Program

The first of these tools are a set of WinWedge configuration files that are designed to make it easy to input weight values directly into any Windows program. The files are named: BalanceCOM1.SW3, BalanceCOM2.SW3, BalanceCOM3.SW3, BalanceCOM4.SW3, BalanceCOM5.SW3, BalanceCOM6.SW3, BalanceCOM7.SW3 and BalanceCOM8.SW3.

To use these files, you will first need to determine which COM port your balance is connected to by looking in the Windows Device Manager in the “Ports COM & LPT” section.

If your balance is connected to COM1, double click on the file named BalanceCOM1.SW3 (likewise, if your balance is connected to COM6, you would double click the file named BalanceCOM6.SW3). This will launch and activate WinWedge on COM port and it will also minimize WinWedge to the system tray to get it out of the way.


From this point on, you should be able to open any Windows program and when you press the function key F9 on your keyboard (or press the Print button on your balance), you should see a weight value pop into whatever program has the input focus wherever the cursor is.

Collecting Balance Readings Using a Ready-Made Excel Workbook

Also included with the Sartorius-Wedge software package is an Excel workbook named SartoBalance.XLSM. SartoBalance.xlsm is a simple balance data collection workbook that will read in data from any Sartorius balance and stack weight values up in a column along with date and time stamps for each reading. The workbook will also chart the data as it is being received as well as perform some basic statistical calculations on the collected data. It can also be used to collect balance readings at a user defined rate (up to 10 times a second).

To use the SartoBalance.XLSM workbook, first make a note of which COM port your balance is connected to and then open the SartoBalance.XLSM file in Excel. If you do not know what COM port your balance is connected to, you should be able to look in the Windows Device Manager under the “Ports COM & LPT” section to determine what COM ports are installed in your system.

Once you know the COM port that the balance is connected to, all you need to do is to open the SartoBalance.XLSM workbook and choose the COM port from the Setup window that is displayed when you open the workbook and then click the OK button in the Setup window. When you do this, WinWedge should automatically launch and activate itself on the selected COM port and you can then use buttons on the left side of the workbook to input data from your balance.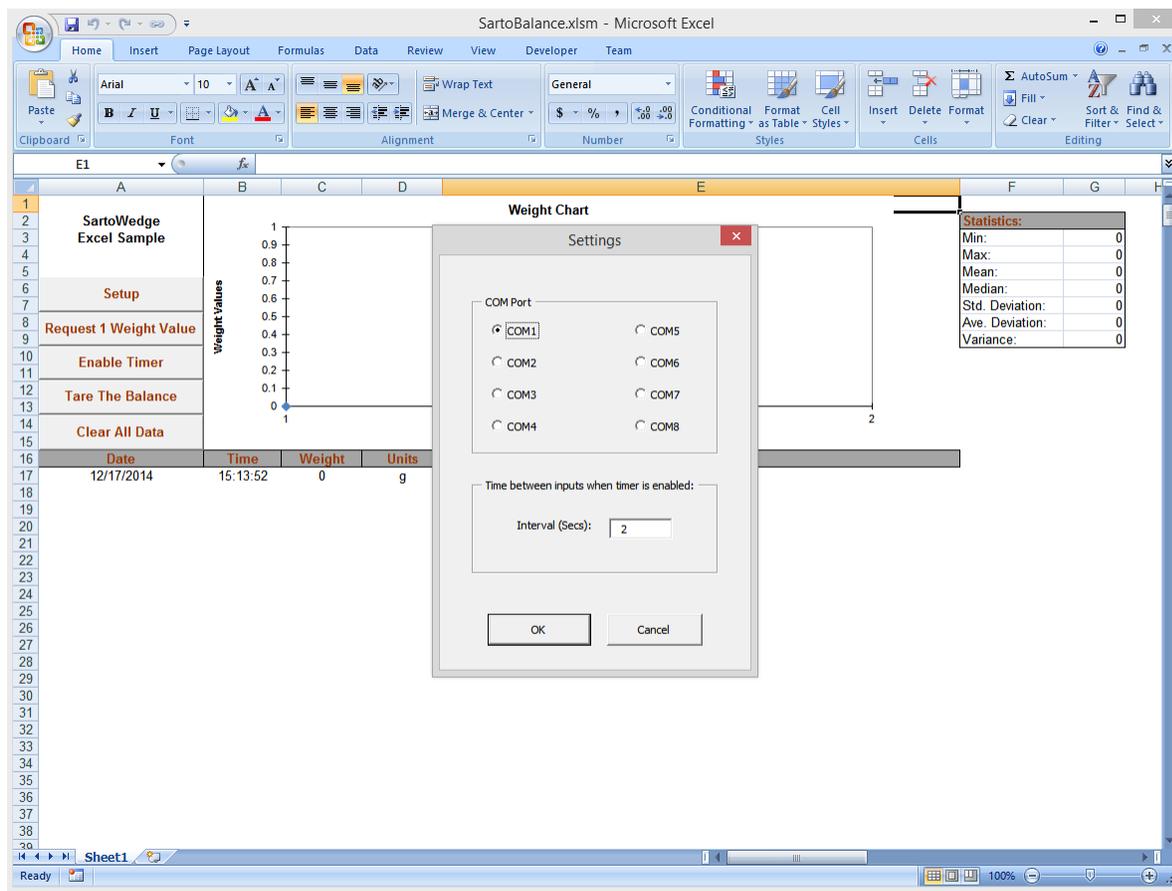


Inputting Balance Readings into Any Excel Workbook


The file named SartoBalance.XLAM (included with the Sartorius-Wedge software) is an Excel file that is designed to be installed as an Add-In to Microsoft Excel so that you can input weight readings from a balance into any open Excel workbook.

To install the SartoBalance.XLAM Add-In, run Microsoft Excel and select: File – Options – Add-Ins and then when the Add-Ins window opens, click the button labeled “Go...” at the bottom of the window next to the “Manage” and “Excel Add-Ins” options.

The following is a screen shot of the Excel Options window: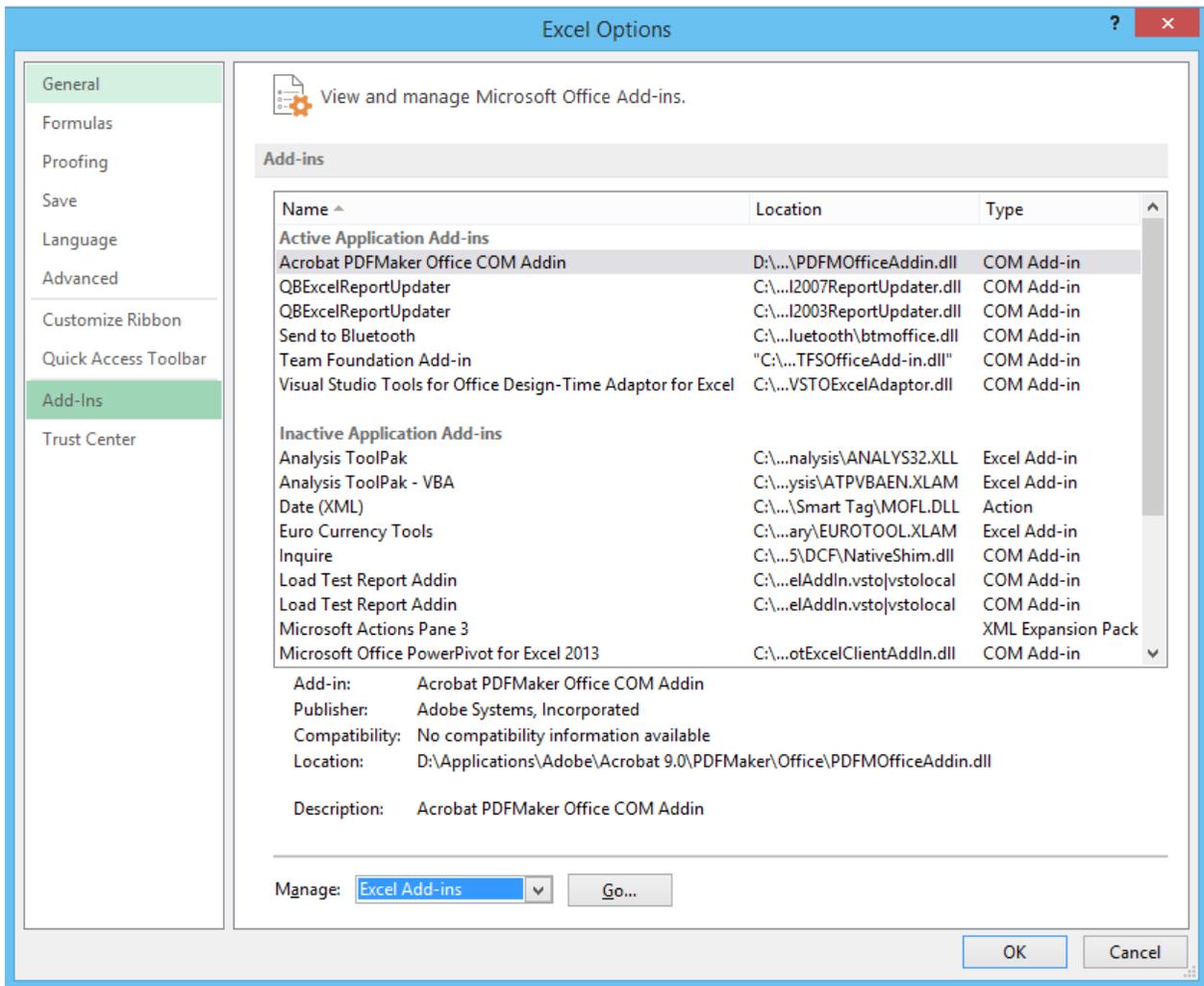

Clicking the “Go...” button will open the following window where you can Browse for Add-Ins.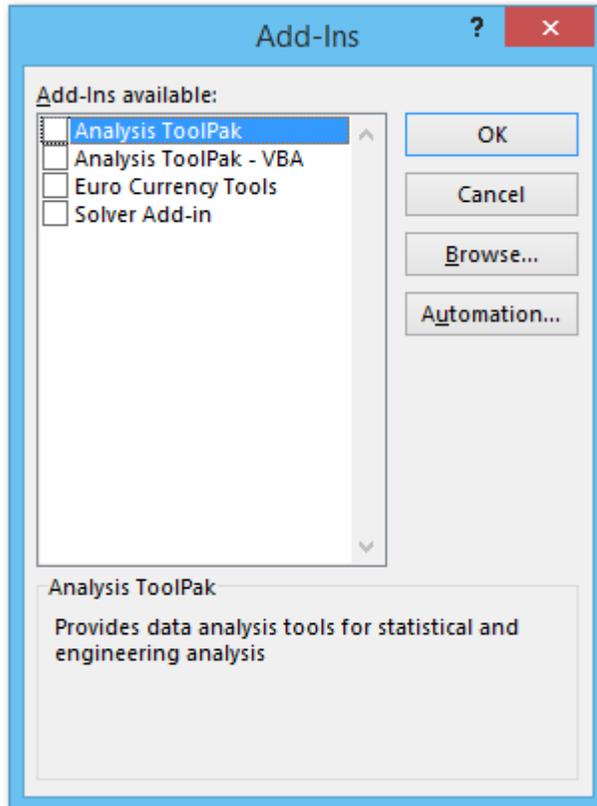

Click the “Browse” button and then navigate to the folder where the Sartorius-Wedge files are installed in your system and select the file named SartoBalance.XLAM. The file should appear in the list of available Add-Ins with the name SartoBalance. Make sure that the checkbox next to the SartoBalance add-in is checked and click the OK button.

From this point on, no matter what Excel workbook is open, there should be an Add-Ins menu in the Excel main toolbar and if you select the Add-Ins menu, you will see a “Sartorius-Wedge” drop down menu with the options “Setup”, “Input 1 Weight”, “Enable Timer” and “Tare the Balance”.

Before you can input weight readings, you need to click the “Setup” menu option and select the COM port that the balance is connected to and click OK. This will launch and activate WinWedge for the COM port that the balance is connected to. Once WinWedge has been launched, you should be able to put the cursor in any cell in whatever Excel workbook is open and when you press function key F9, you should see a weight reading appear in the active cell and the cursor should move down one cell.

Even if you close Excel, when you re-open it, the Add-In will still be available in all workbooks. To remove the add-in, go through the exact same steps as above and remove the check mark next to the SartoBalance Add-In in the list of available Add-Ins.

Note, you cannot use the SartoBalance.XLSM file at the same time as the SartoBalance.XLAM Add-In. If you want to use the SartoBalance.XLSM workbook, you must first remove the SartoBalance Add-In by removing the check mark next to the SartoBalance Add-In in the list of available Add-Ins.

Connecting a balance to an Ethernet network

To connect your balance to an Ethernet network, you will need to understand how TCP/IP network devices work and how they communicate with each other.

Every device connected to an Ethernet network has a unique “IP Address” that is similar to the main telephone number for a company. Just like a telephone call, one side of the connection (called the “Client”) needs to know the IP Address of the device that it wants to connect to and it also needs to know the specific “Port” number to connect to. The Port number is similar to a telephone extension number. The device that is waiting for calls to come in is called the “Server” and it must be configured to “listen” for connections on a specific TCP/IP port number and accept connection requests when they come in on that port number. Some port numbers are reserved for specific jobs like serving up web pages or sending or receiving email however most port numbers can also be configured to just send or receive raw data similar to a RS232 serial port. Port numbers can range from 1 to 65535.

The IP Address that a device uses can either be assigned automatically by the router that you connect the balance to or you can also hard code the IP Address into the device. An IP address that is assigned by a router is called a “Dynamic IP Address” and a hard coded IP Address is called a “Static IP Address”.


In the case of a Sartorius balance, you can configure it to either use a Dynamic or a Static IP address. The problem with Dynamic IP Addresses is that you are not guaranteed to get the same IP Address assigned to the device each time you power it on. Because of this, you are typically better off setting a static IP Address for the balance. The Port Number that is used for the connecting to the balance is also configurable in the balance network settings menu.

For the physical connection between the balance and your network, you should use a standard Ethernet cable and connect it to either a router or a switch in your local area network.


Example - Configuring the Network Settings in a Cubis Balance

To configure a Cubis balance network settings, you would select the “Menu” button on the balance display and then select “Device Parameters”, “Configure Ports” and “Network Settings (Ethernet)”. This will display the following screen where you can set up the balance to use a fixed IP Address and also enter the IP Address that you want the balance to use. You may need to consult with your network administrator to select an appropriate IP Address that fits with your particular network.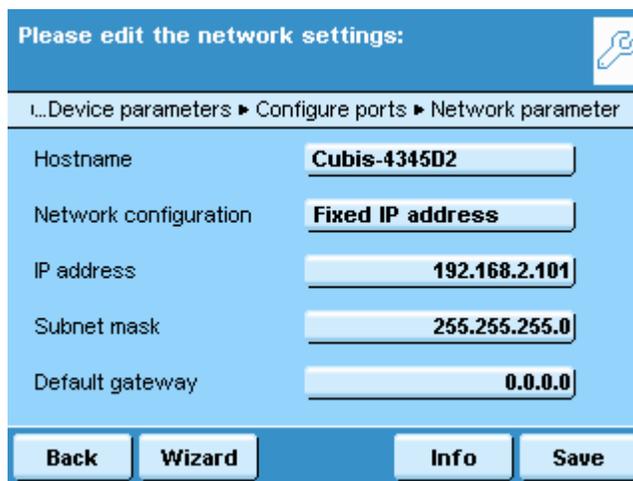

Please edit the network settings:	
...Device parameters > Configure ports > Network parameter	
Hostname	Cubis-4345D2
Network configuration	Fixed IP address
IP address	192.168.2.101
Subnet mask	255.255.255.0
Default gateway	0.0.0.0
Back Wizard Info Save	

Once the IP Address is set, you can click the “Save” button to save your settings and then click the “Back” button to return to the previous menu where you can select the option “Network port setup”. When you do so, the following screen will appear: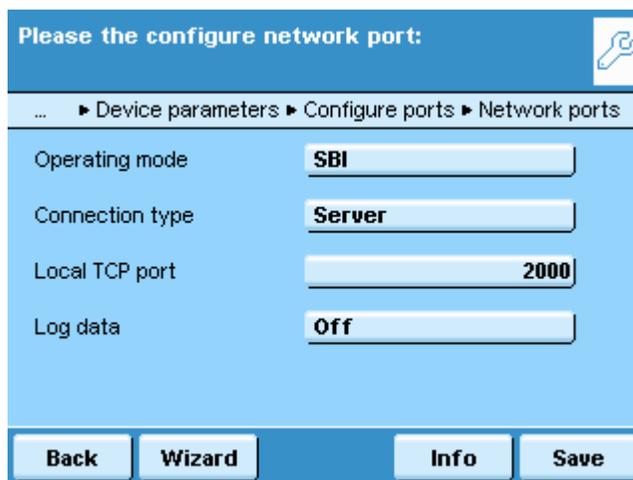

Please the configure network port:	
... > Device parameters > Configure ports > Network ports	
Operating mode	SBI
Connection type	Server
Local TCP port	2000
Log data	Off
Back Wizard Info Save	

In the above screen, you should set the “Operating Mode” setting to “SBI (Sartorius Standard Protocol)” and set the “Connection Type” to “Server” and then enter the “Local TCP Port” that you would like to use. You can enter any number that you like below 65535 however we recommend that you select a number above 1000. In this example, we have set the Local TCP Port to the value 2000.


When finished, click the “Save” button to save your settings and then click the “Back” button until you return to the main balance display screen where the weight appears.

At this point the balance is configured as a TCP/IP Server with a static IP Address and it is waiting for a TCP/IP Client to connect to it on port number 2000. The next step is to configure the TCPWedge software (included with Sartorius-Wedge) to use these connection settings.


Configuring TCPWedge to Connect to the Balance

The TCPWedge program included with the Sartorius-Wedge package is a special version of the WinWedge software that is designed for connecting to network devices instead of RS232 Serial devices.

When you run TCPWedge.exe, the following main menu will appear: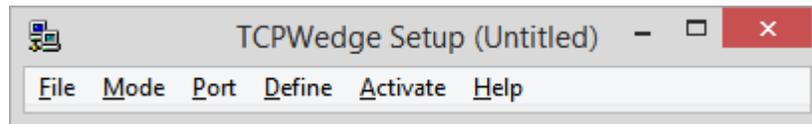

Typically, the first step in configuring TCPWedge is to choose how you want the balance data transferred to another program and optionally what program to transfer it to. This is done by selecting the "Mode" menu. The simplest way to transfer data to another program is by sending the data as keystrokes. From the Mode menu, select the option "Send Keystrokes To" and then in the window that appears, clear out the textboxes labeled "Application Title Bar Text" and "Command Line". Clearing out these two items will cause TCPWedge to send your weight values to whatever application has the input focus when the weight data is sent from the scale. The window should appear as shown below. Click the OK button to return to the TCPWedge main window.

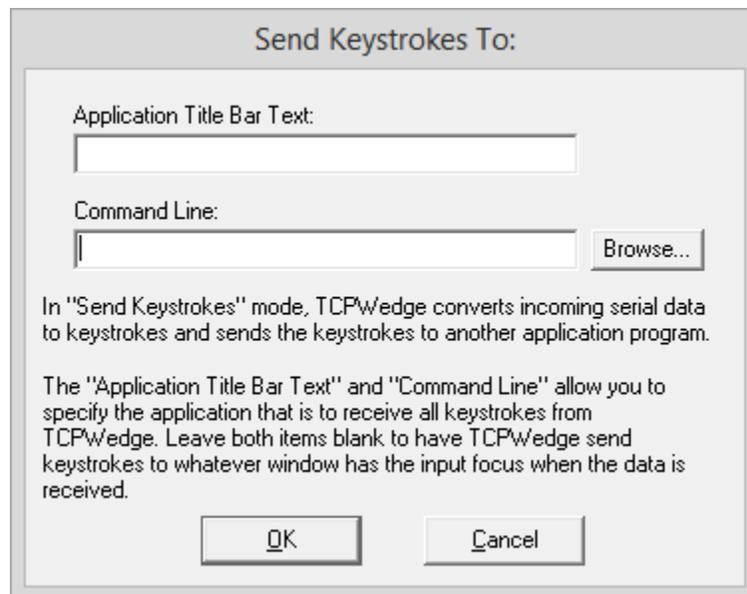The next step is to select "Port" and "Settings" from the main menu. This will open the following window where you can enter the IP Address of your balance and the TCP/IP port number that you configured it to use. In this

example, we would enter the IP Address 192.168.2.101 and the TCP/IP Port number 2000 as shown below. When finished, click the OK button to return to the main menu.

The screenshot shows a dialog box titled "TCPWedge Port Settings". It contains two radio buttons: "This PC will act as TCP/IP Client" (which is selected) and "This PC will act as TCP/IP Server". Below these are two text input fields: "Remote Host IP Address" with the value "192.168.2.101" and "Remote Port" with the value "2000". At the bottom are "OK" and "Cancel" buttons.


The Sartorius balance transmits weights as a 22 byte string containing the weight value and one or two characters indicating the units (i.e. g for grams, oz for ounces, etc.). In most cases you will only want the numeric weight value sent to another application followed by an Enter or a Tab keystroke. To configure TCPWedge to work this way, select "Define" - "Input Data Record Structure" from the main menu and then click the Continue button in each window that appears until you get to the "Input Record Definition Editor - Send Keystrokes Mode" window. Change the "Input Filter" option to "Numeric Data Only" and then click your mouse in the textbox labeled "Field Postamble Keystrokes" and then click the button labeled "Keystroke List". In the list of keystrokes that appears, select the "Enter" keystroke and click OK. The window should appear as below. Click OK to return to the main menu.

The screenshot shows a dialog box titled "Input Record Definition Editor - Send Keystrokes Mode". It features a "Record Preamble Keystrokes" text box at the top. Below is a table with four columns: "Field", "Input Filter", "Math Expression", and "Field Postamble Keystrokes". The "Field" column contains the number "1". The "Input Filter" column has a dropdown menu set to "Numeric Data Only". The "Field Postamble Keystrokes" column has a text box containing "{ENTER}". At the bottom are "OK", "Keystroke List...", and "Cancel" buttons.

Field	Input Filter	Math Expression	Field Postamble Keystrokes
1	Numeric Data Only		{ENTER}


The Sartorius balance will transmit a weight value when you press the “Print” button on the balance and you can also trigger the balance to transmit a weight by sending the Escape character (ASCII 27) followed by an upper case P. WinWedge and TCPWedge both have the ability to define hotkeys that you can use to send character strings out the port to the balance. This means that you can configure a hotkey on your keyboard to trigger the balance to send a weight reading when you press the hotkey.

To configure a hotkey in TCPWedge, select Define – Hotkeys and Hotkey Actions from the main menu. In the window that appears, select the option “Transmit String” from the list of Hot Key Actions and then put the cursor in the textbox labeled Hot Key Keystroke and press function key F9 on your keyboard. Next, put the cursor in the textbox labeled “Output String” and then click the button labeled “ASCII Chart” and select the ASCII 27 character (Esc) from the list of characters and click the OK button. Next, type an upper case P on your keyboard. The window should appear as it does below. Click the OK button to return to the main menu.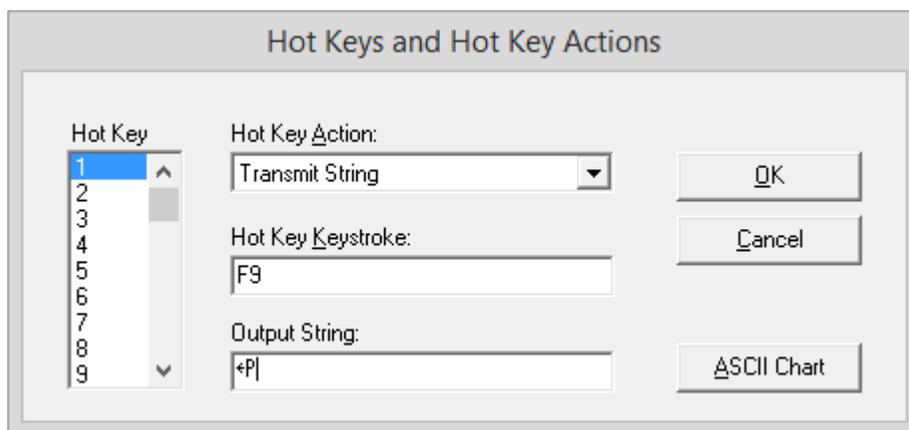


At this point, TCPWedge is fully configured to input data from your balance into any Windows program. Select “File” – “Save” and save your settings to a TCPWedge configuration file so that you do not have to go through the above steps to configure TCPWedge again. If you save your configuration file to your Windows desktop, you can launch and activate TCPWedge just by double clicking the configuration file icon on your desktop.

Once your settings are saved, the last step is to activate and test TCPWedge. This is done by selecting “Activate” – “Normal Mode” from the main menu.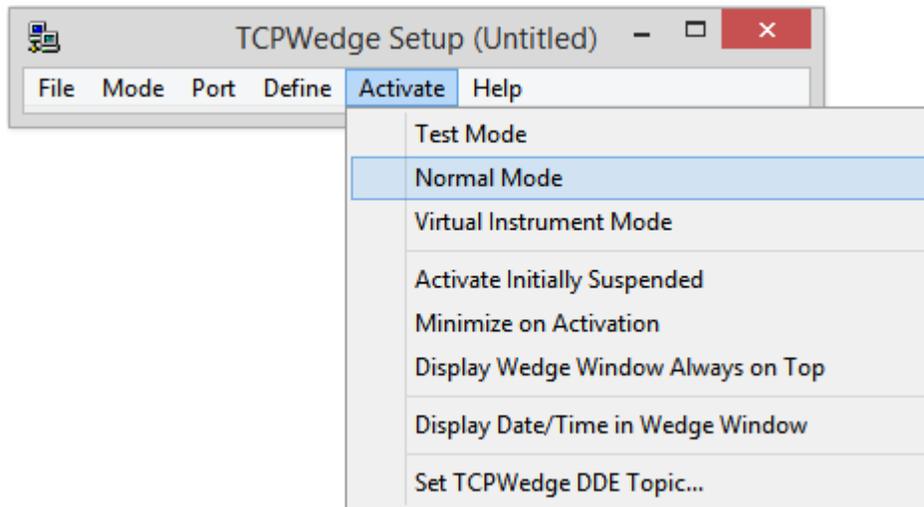

When you do this, you should see the window shown below.

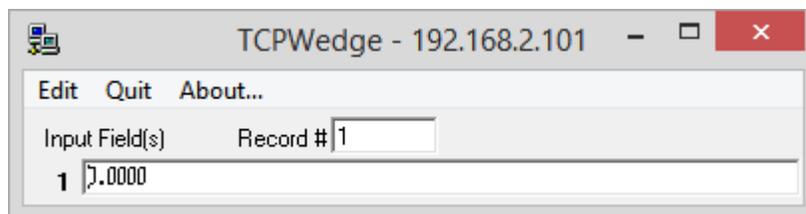At this point, you should be able to open any Windows application that you like including Excel, Word, Notepad, Access, etc. and when you press function key F9 on your keyboard, a weight value should appear wherever the cursor is in the other application. You can also press the Print button on your balance to transmit a weight reading into the application.

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